



## PRODUCT INSERT

# Nutri-Fly™

## Instant Formulation

Powdered Drosophila Media by Genesee Scientific

**Description:** Nutri-Fly Instant Formulation is made to accommodate the requirement of labs without the ability to cook traditional Drosophila media or for labs making very small amounts (<1L) of Drosophila media at a time.

**Directions:** Shake the bag of media prior to measuring to ensure a more even distribution of ingredients. Add equal parts Nutri-Fly Instant Formulation and water to your vial or bottle. Mix to ensure all dry media comes in contact with water. Let medium sit for 5 minutes prior to adding flies.

**Ingredients:** potato flakes, dextrose, brewer's yeast, tegosept

Cat#	Makes
66-117	~1L
66-118	~4L

### **Helpful Hint**

The proper ratio of water to Nutri-Fly Instant Formulation may vary depending on your particular environment or the size of your culture vessel. Generally, you should use less water than the volume of instant media for larger vessels (<6oz). If this is your first time using Nutri-Fly Instant Formulation, test a small amount first to find the ratio of media to water that works best for you.